

SEDEX Low Temperature Evaporative Light Scattering Detector

SAMPLE: POLYETHOXYLATED ALCOHOLS

HPLC

SFC

POLYETHOXYLATED ALCOHOLS MIXTURE



Column Lichrospher 100 DIOL 125 x 4 mm 5 μ m
ELSD 45°C

HPLC

A. hexane
B. CHCl₃ + Isopropanol 2%
Flow rate: 1 ml/min

SFC

CO₂/MeOH 87/13 w/w BC20.BC40
CO₂/MeOH 93/7 w/w BC30

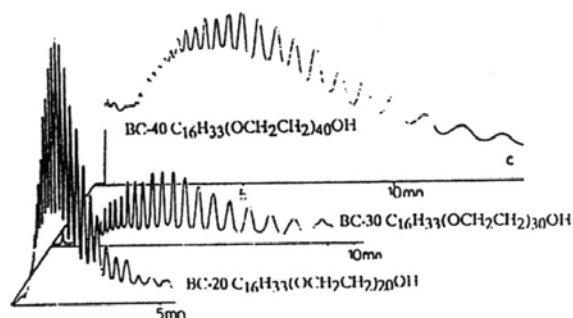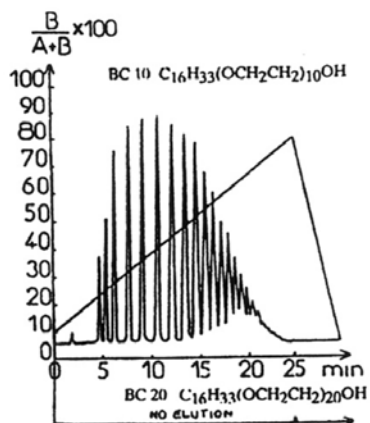

Analyzed by: S. Brossard, M. Lafosse and M. Dreux [J. Chromatogr. 591 (1992) 149-157]